SPECTROLINE® FC-100 Series Fan-Cooled, Long Wave UV Lamps

Ultraviolet radiation emitted from this product. Avoid exposure. ALWAYS WEAR PROTECTIVE CLOTHING. EXPOSURE MAY CAUSE PREMATURE AGING OF THE SKIN AND CANCER. ALWAYS WEAR PROTECTIVE EYEWEAR; FAILURE TO DO SO MAY RESULT IN SEVERE BURNS OR LONG TERM INJURY TO EYE. Never look directly into the lamp. Exposure can cause eye and skin allergy and allergic reactions. Medications or cosmetics may increase your sensitivity to ultraviolet radiation. Consult physician before operating this product if you are using medications or have a history of skin problems or believe yourself especially sensitive to sunlight. FOR PROFESSIONAL EXAMINATION USE ONLY

IMPORTANT SAFETY INSTRUCTIONS

1. Read all instructions.
2. Care must be taken, as burns can occur from touching hot parts.
3. Do not operate the lamp with a damaged cord or if the lamp has been damaged.
4. **Do not look directly into the UV light. Wear UV-absorbing protective eyewear (such as the UVS-30 glasses provided) when operating this or any UV lamp. Never operate the lamp with the UV filter lens removed.**
5. Never operate this equipment in any manner not specified in these instructions because the protection provided by the equipment may be impaired.
6. Adequate ventilation should be provided when working on operating combustion engines.
7. This unit is not recommended for high-dust environments.

GENERAL INFORMATION

1. Install the UV bulb and the Bulb-Saver™ (packed separately).
2. Connect the plug to a power outlet that complies with the electrical requirements specified on the lamp's label.
3. The **120V** and **100V** lamps require just **one three-position switch** (on the ballast) to operate. The first position (*lowest*) is OFF. The second position (*up*) turns on the fan, the third position (*up*) turns on the lamp.
4. The **230V** and **240V** lamps require **two switches** to operate. Turn ON the ballast switch. This powers both the fan and the lamp head. **The unit will not operate unless the ballast/fan switch is turned ON first.** Turn ON the switch on the lamp handle. Allow approximately 3 to 5 minutes for the lamp to reach full lighting intensity. If the lamp fails to light, check that the bulb is completely screwed into the socket and that both the switches are turned ON.
5. When the lamp is used intermittently and frequently, do not turn it off until the inspection is complete. Any interruption in the power supply, such as a drop in voltage or changing power outlets, will cause the bulb to go out. The bulb must then cool for a minimum of 5 minutes before restarting. Continuous operation will also extend the bulb life.
6. Turn OFF the switch on the lamp handle, **but leave the ballast/fan switch ON for 4 to 5 minutes** to reduce heat buildup in the lamp head. Turn off all power before storing the lamp.
7. **CAUTION: The lamp housing is fan-cooled, but the filter lens will become hot during operation. The lamp stand is attached to the unit for increased protection. Do not remove it during operation. Do not stand the lamp on furniture or papers while in operation.**

OPERATOR SERVICING

Do not perform any servicing that is not covered in these instructions. Return the lamp to the factory for maintenance not covered in this section.

HOUSING AND FILTER MAINTENANCE

- Always unplug the lamp from the power outlet before performing any maintenance or servicing. Allow the UV lamp to cool before handling.
- Immediately clean spills from the housing with water and a mild detergent and wipe dry. Periodically clean the filter with a glass cleaner and a soft cloth.

BULB/FILTER INSTALLATION OR REPLACEMENT

Use only a Spectrolines® 100S/M UV bulb. Using any other bulb may damage the unit and will void the warranty.

1. Remove the lamp stand/protector.
2. Detach the silicone-rubber filter holder by peeling it from around the circumference of the lamp housing. The filter will remain inside the filter holder. (If replacing the filter, press it firmly but gently inside the filter holder.) Set the filter holder aside.
3. Place the UV bulb through the Bulb-Saver™/collar in the lamp housing. Turn the bulb *clockwise* gently but firmly into the socket with the fingertips. (If removing an existing bulb, turn it *counterclockwise*.) **Do not operate the lamp without the Bulb-Saver.**
4. Reattach the filter holder by pressing it with the heel of your hand around the circumference of the lamp housing. Be sure the groove in the filter holder seats over the rim of the housing.
5. Reattach the lamp stand/protector.

UV INTENSITY VERIFICATION

To ensure that your lamp is operating at the required UV intensity, it should be checked periodically. The AccuMAX™ XRP-3000 radiometer/photometer kit uses a dual-wavelength UV-A/VIS sensor detector to accurately measure ultraviolet irradiance or visible light. The AccuMAX XR-1000 digital readout unit with the XS-365 UV sensor or the Spectrolines DM-365XA digital readout unit are also recommended for accurate UV-A measurement. The meters are specially designed to measure the UV irradiance from 320–400nm with a peak at 365nm.

For more information about these units, please contact the Customer Service Department at Spectronics Corporation. In the U.S. and Canada, call toll-free 1 800-274-8888. Outside the U.S., call 516-333-4840.

ELECTRICAL SPECIFICATIONS

Model	Filter	120V/60Hz	230V/50Hz	240V/50Hz	100V/50-60Hz
FC-100	2F958	1.5 Amps	—	—	—
FC-100D	2F350	1.5 Amps	—	—	—
FC-100/F	2F958	—	0.9 Amps	—	—
FC-100D/F	2F350	—	0.9 Amps	—	—
FC-100/FA	2F958	—	—	0.95 Amps	—
FC-100D/FA	2F350	—	—	0.95 Amps	—
FC-100/FB	2F958	—	0.9 Amps	—	—
FC-100D/FB	2F350	—	0.9 Amps	—	—
FC-100/J	2F958	—	—	—	2.7 Amps
FC-100D/J	2F350	—	—	—	2.7 Amps

All bulbs are 100S/M, 100W Medium Base Spot

The FC-100D Series lamps are equipped with a diffusing filter that produces a "diffused spot" beam of UV-A light.

ENVIRONMENTAL CONDITIONS

The FC-100 Series UV lamps are designed to be safe under the following conditions:

- Indoor use;
- Altitude up to 2,000 m (6,562 ft.);
- Temperature 5°C to 40°C (41°F to 104°F);
- Maximum relative humidity 80% for temperatures up to 31°C (88°F) decreasing linearly to 50% relative humidity at 40°C (104°F);
- Mains supply voltage fluctuations not to exceed ±10% of the nominal voltage;
- Installation Category II;
- Pollution Degree 2.

WARRANTY

The warranty policy for the FC-100 Series UV lamps is provided on the Certificate of Limited Warranty enclosed separately with each unit.

NOTE: For assistance of any kind, contact the Customer Service Department at Spectronics Corporation.

Item	Part No.	Description
1	FP-175	Filter Protector/Lamp Stand
2	123604	Filter Holder
3	2F958	Filter Lens
	2F350	Filter Lens, Diffusing
4	100S/M	Bulb, Spot
5	106862	Bulb-Saver™
6	EC-0415	Socket, Medium Base
7	124736	Lamp Housing
8	107394	Socket Plate
9	122833	Adapter, Fan
10	123026	Fan, 120V, 100V
	123027	Fan, 230V, 240V
11	122832	Fan Cover
12	100310	Pistol Grip
13	118677	Handle, Ballast
14	108291	Ballast Assembly (120V)
	105049	Ballast Assembly (230V, 240V)
	117479	Ballast Assembly (100V)
15	108388	Power Switch (120V, 100V)
	100253	Power Switch (230V, 240V)
16	119797	Cord, Primary, 8 ft
	119798	Cord, Primary, 20 ft
17	WC-0130	Cord, Secondary
18	101145	Power Switch, Handle (230V, 240V)

One switch is required on 120V and 100V units. This switch is located on the ballast assembly. Each of the primary cords is fitted with the appropriate plug for the country of shipment: Domestic, 120V; European, 230V; British, 230V; Australian, 240V or Japanese, 100V. For additional information, contact Spectronics Corporation.

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HINWEIS: Wenden Sie sich zwecks etwaiger Hilfestellung bitte an die Kundendienstabteilung der Spectronics Corporation.
PRINTED IN U.S.A.